


TGA Balance Purge Switch

For Simple Gas Changeover for TGA Without
Using Any Tools or Undoing Gas Connections

www.lab-ware.co.uk



Changing gases to any analyser usually involves taking out some tools and opening connections to change to a different gas. This can involve shutting down the gas system to avoid wasting gas, or may cause leaks when reconnected. The RSL gas switch removes the need to change any connections, and instead a simple changeover valve switches the gas to the new configuration. The existing second gas outlet is not affected by the switch, although different configurations can be provided as a customisation, usually at a small additional cost.

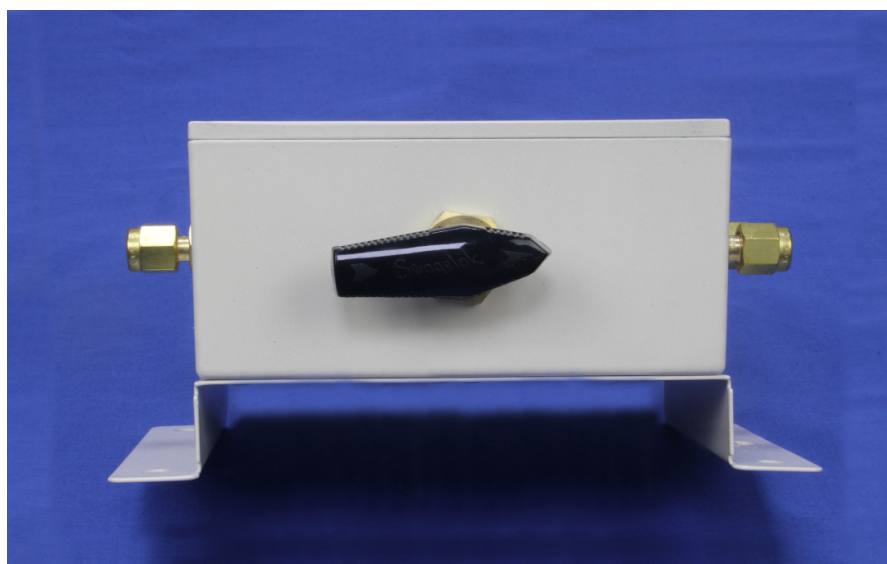

Ruston Services Ltd, 9 Broomfield Gardens, Shandon, Helensburgh, Argyll & Bute, G84 8HR, United Kingdom

Telephone +44 (0) 7980 222573

Email info@lab-ware.co.uk

Website www.lab-ware.co.uk

TGA Balance Purge Switch

www.lab-ware.co.uk

Specifications

Case	All-Metal case with four 5mm wall mounting holes at 76mm x 144mm centres.
Inlet Connections	Brass Swagelok™ connection for 1/8 inch pipe connections
Outlet Connections	Brass Swagelok™ connection for 1/8 inch pipe connections
Internal Pipework	All in metal tubing to avoid gases diffusing from outside the pipes
Dimensions	120mm x 165mm x 78mm (W x H x D)
Shipping Weight	0.70 kg
Configuration Options	Different configurations of the inlet and outlet connections are available - please inquire.

Ordering Information

Quotation Requests	Please email your requirements to sales@rustons.org.uk , stating what you would like to order Quotations are normally valid for 60 days.
Ordering	Please send your official order number to the address below or by email to sales@rustons.org.uk If your organisation is VAT registered, please ensure that your VAT number is clearly visible on your order.
Delivery	Delivery is normally 4 - 6 weeks from receipt of order. We may be able to provide faster delivery in some circumstances



Ruston Services Ltd, 9 Broomfield Gardens, Shandon, Helensburgh, Argyll & Bute, G84 8HR, United Kingdom

Telephone +44 (0) 7980 222573

Email info@lab-ware.co.uk

Website www.lab-ware.co.uk